

Dino-Lite AMK4012-C200

Dino-Lite AMK4012-C200 at a glance



AMK-4012

In the box



C200



TV Connector










Battery

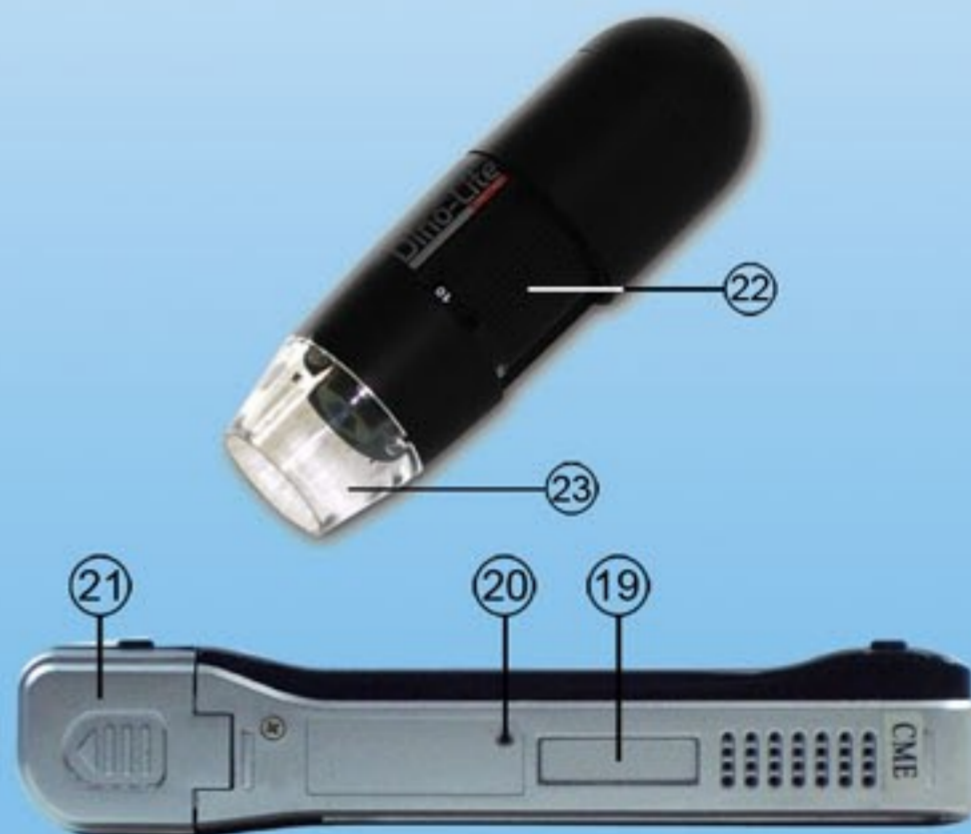

Battery Charger

Function Keys



1. Power Switch	9. Volume Control 	17. A/V Recording  (REC)
2. Speakers	10. Remote Control Receiving Interface	18. Key Lock
3. Charging Status Indicator	11. SD Card Slot	19. Expanding Slot Connector
4. Recording Status Indicator	12. USB Slot	20. Reset Key
5. Setup 	13. Battery Charging Slot	21. Battery Cover
6. Snap 	14. Microphone	22. Focus & Magnification Adjustment Knob
7. Stop 	15. AV-Input	23. LED Lighting & Cover
8. 5-way Navigation Keys 	16. AV-Output / Earphone jack	

Dino-Lite AMK4012-C200

Charging the Battery of C200

For the first time user, please charge the C200 battery for more than 8 hours by connecting the C200 to the power adapter and then connect to an power outlet. You can also charge the C200 with an USB cable (sold separately).

Notice:

Even when the adapter is connected, the C200 will not start when the battery is discharged.

To extend the battery life, please unplug the adapter immediately when done charging the battery.


Starting up the AMK4012-C200

1. Connect Dino-Lite AMK4012 to AV-Input of C200 before powering on the C200.
2. Make sure the SD card is inserted in the SD card slot of C200.
3. Turn on the C200 by pressing the power switch. The LED of AMK4012 will turn on, and the following screen shown below will be displayed initially for approximately 2 seconds. After that, the screen will then display the real-time video transmitted from the AMK4012.


Taking Picture



Taking a picture can be done by simply pressing .
The picture will be stored in the SD card.

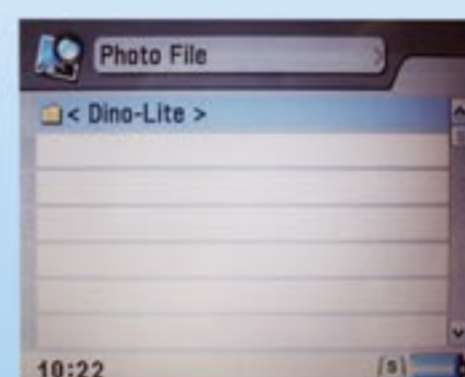

To review the picture, press one time to enter the Photo File. The <Dino-Lite> folder can then be found.

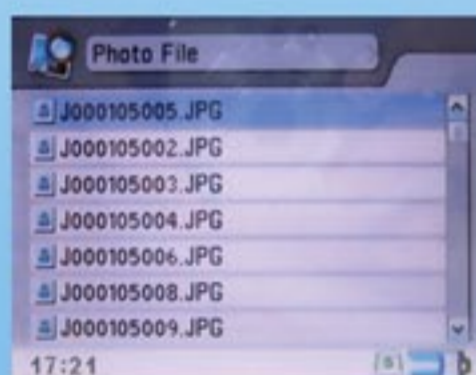


Access the <Dino-Lite> folder by pressing , and a list of picture files will be shown as above.



Scroll through the Photo File list by using to the file that you want to select, then press and the picture will be displayed.




The picture can be rotated 90-degrees by pressing .

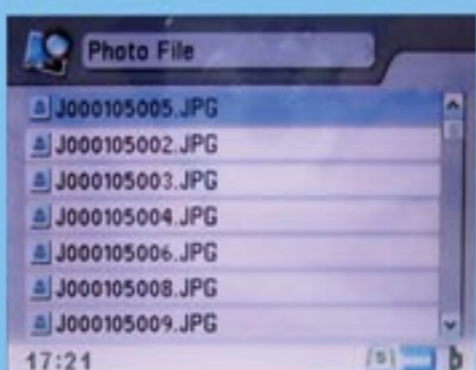

Return to the Photo File list or the preview screen by pressing .



Delete the picture in the Photo File list by pressing and the following screen will be displayed.

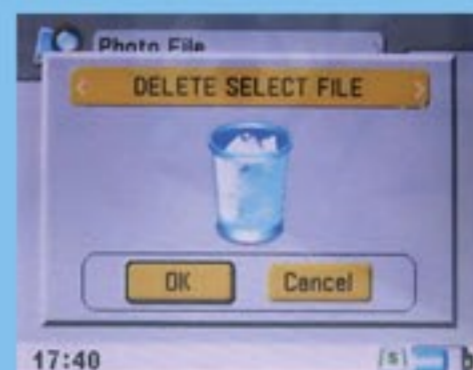


Delete a single picture or all pictures by selecting One or All in the delete option.

Taking Video




Taking a video can be done by simply pressing the Record button  (REC). To pause the recording, press the Record button  (REC) again. Stop or return from the recording by pressing . While recording a video, audio is also recorded with its built-in microphone. The recorded video will be stored in the SD card.


To review the video, press  **two times** to enter the Video File window. The <Dino-Lite> folder can then be found.




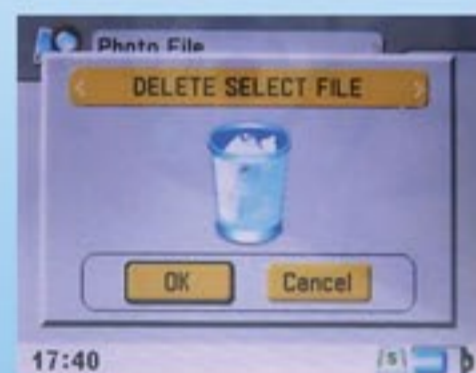
Access the <Dino-Lite> folder by pressing , and a list of video files will be shown as above.

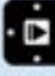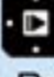

Scroll through the Video File list by using [ : up/down] to the file desired then press  and the video will be displayed. Return to the Video File list or the preview screen by pressing .

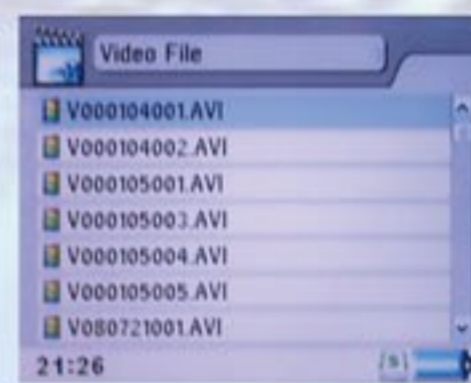


Deleting a video in the Video File list can be done by pressing  and the following screen will be displayed.




Delete a single video or all videos by selecting One or All in the delete option.

Setup

The setup screen provides functions including: time setting, TV format, display, language settings, format the SD card, backup from SD to harddisk (if a harddisk is available), firmware upgrade etc.

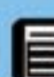
By pressing  **three times** in the preview mode, a setup screen will then be displayed as following.

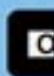

Use  to select the Setup menu and change it's setting. Exit Setup by pressing  simply wait approximately 3 seconds and the C200 will automatically exit the Setup menu.

Connecting the AMK4012

Connect to AV

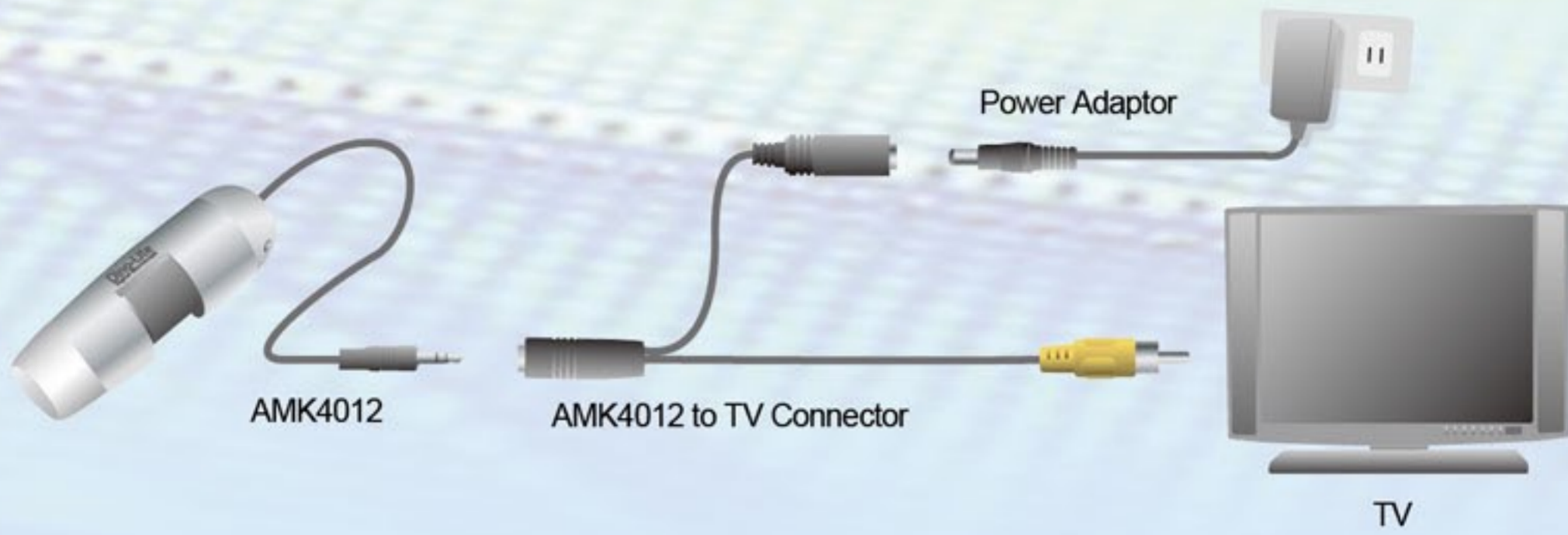

Connect to C200



SD Compatibility

- SD card up to 16GB class 6
- Also Eye-Fi™ SD card compatible (SD card not included)

